

## Operation of Chameleon UltraII IR laser

**Note:** The water recirculator is **ALWAYS ON**. The temperature should read **20 C**  
Refer to diagram below for Power supply/Control box switches and buttons

1. Confirm that the AC ON light (1) is lit.
2. Turn key (2) from STANDBY to ON.
3. The Display (4) will provide status information: the last wavelength, power and will indicate OK on bottom of display when the laser is modelocked at the indicated wavelength.
4. To change wavelength, turn the POWER ADJUST knob (11) to desired wavelength and then push the MENU SELECT button (6)
5. Once the laser is ready, open the SHUTTER (8)
6. The laser will operate at full power, there is no way to adjust the output, other than by wavelength.

***BE CAREFUL TO NOT PUT EYES AT BEAM LEVEL! IF YOU ARE GOING TO DIVERT THE BEAM, WEAR APPROPRIATE EYE SAFETY GEAR (on shelf).***

To Turn **OFF** - Close Shutter (8) and turn key (2) to STANDBY.

**If you have trouble, please seek assistance.**

